2000005144

UTTD_052016

IKA®

IKA® ULTRA -TURRAX® Tube Drive

Betriebsanleitung	DE	4
Ursprungssprache		
Operating instructions	EN	11
Mode d'emploi	FR	18
Руководство пользователя	RU	25
Indicaciones de seguridad	ES	32
Veiligheidsinstructies	NL	33
Norme di sicurezza	IT	34
Säkerhetsanvisningar	SV	35
Sikkerhedshenvisninger	DA	36
Sikkerhetsanvisninger	NO	37
Turvallisuusohjeet	FI	38
Instruções de serviço	PT	39
Wskazówki bezpieczeństwa	PL	40
Bezpečnostní pokyny	CS	41
Biztonsági utasítások	HU	42
Varnostna opozorila	SL	43
Bezpečnostné pokyny	SK	44
Ohutusjuhised	ET	45
Drošības norādes	LV	46
Nurodymai dėl saugumo	LT	47
Указания за безопасност	BG	48
Instrucțiuni de securitate	RO	49
Υποδείξεις ασφαλείας	EL	50

IKA-Melco, Germany
Reg. No. 004343

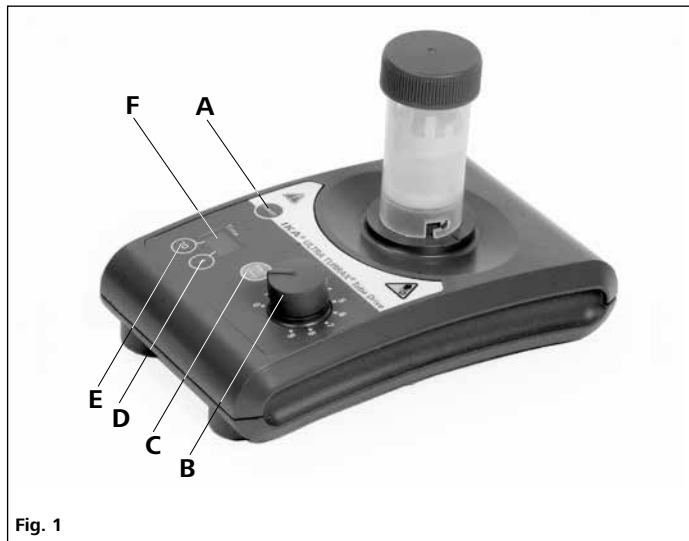
Fig. 1

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Contents

	Page
Declaration of conformity	11
Explication of warning symbols	11
Safety instructions	12
Unpacking	13
Correct use	13
Useful facts	13
Commissioning	13
Error message	15
Accessories	15
Maintenance	16
Warranty	16
Technical data	17

Declaration of conformity

We declare under our sole responsibility that this product corresponds to the regulations 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the standards or standardized documents EN 61010-1, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

Explanation of symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information can result in damage to the device or to system components.

Risk of hand injury.

Safety instructions

For your protection

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- The device must be supervised at all times when in operation.

- **Caution!** For research purposes only! Not suitable for diagnostic use (in accordance with the IVD Directive).

DANGER

The device must only be operated with a tube attached; with no tube there is a risk of injury from the rotating locating pin.

- Hearing protection must be worn when working with **BMT-20** or **BMT-50** tubes.

DANGER

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed.

Otherwise there is a risk of:

- splashing liquids
 - projectile parts
 - body parts, hair, clothing and jewellery getting caught.
- Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
 - The feet of the device must be clean and undamaged.
 - The device is not suitable for manual operation.
 - The device may heat up when in use.
 - Check the device and accessories beforehand for damage each time when you use them. Do not use damaged components.

- Ensure that the cover on the tube is screwed on tightly.
- Ensure that the tube is firmly attached to the bayonet lock connector on the drive unit prior to operating the device.
- The tube must only be attached and removed while the motor is stationary.

WARNING

The **IKA**® tubes must always be closed when the device is in operation. Switch off the device immediately if any material leaks from the tube. Clean the device.

- Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation, leading to pressurization of the container: risk of material spraying, protective equipment must be worn.
- The temperature of the material must not exceed 40 °C.
- Only use tubes approved by **IKA**®.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- Beware of the risk of flammable materials.
- **Do not** operate the device in explosive atmospheres, with hazardous substances or under water.
- The device doesn't start up again automatically following a cut in the power supply.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.

For protection of the device

- The voltage stated on the name plate must correspond to the mains voltage.
- The device must only be operated with the original plug-in power supply unit.
- Protect the device and accessories from bumps and impacts.
- The device may only be opened by experts.

Unpacking

- **Unpack:**
 - Please unpack the device carefully
 - In the case of any damage a fact report must be set immediately (for post, rail or forwarder)
- **Delivery scope:**
 - **Tube Drive**
 - Power supply unit
 - Removal hook
 - Operating instructions
 - A warranty card.

Correct use

• Use

	Application	Volume	Mode
▶ with Tube ST-20...	Mixing	2 - 15 ml	Seconds and minutes mode
▶ with Tube ST-50...	Stirring	15 - 50 ml	Seconds and minutes mode
▶ with Tube DT-20...	Dispensing	5 - 15 ml	Seconds and minutes mode
▶ with Tube DT-50...	Stirring	15 - 50 ml	Seconds and minutes mode
▶ with Tube BMT-20...	Crushing	2 - 15 ml	Seconds and minutes mode
▶ with Tube BMT-50...	Mixing wet/dry	15 - 50 ml	Seconds and minutes mode

• Range of use (indoor use only)

- Laboratories - Schools
- Pharmacies - Universities

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network which also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the device is operated with accessories that are not supplied or recommended by the manufacturer
- If the device is operated improperly contrary to the manufacturer's specifications
- If the device is operated improperly contrary to the manufacturer's specifications.

Useful facts

Changes in viscosity and volume caused by processing of dispersion media may cause small fluctuations in the speed of the device.


Commissioning

Pay attention to the ambient conditions listed in the "Technical Data".

Position the rotating knob (B) at the left stop before starting up.

If these conditions are met, the unit is ready for service when the mains plug has been plugged in.

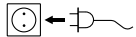
Switching on:

Press "Power" button (A).

Previous mode and time settings will be displayed e.g. :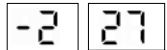


(Factory setting: -2)

Setting Timer (Changing of mode)

Switch off the device.

Minutes-mode


Press and hold "Start/STOP" button (C), and also press the "Power" button (A).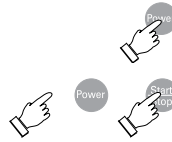

Minutes-mode display:

The time that was previously set will be stored, e.g.:

Note: max. 29 min., speed range upto 4000 rpm.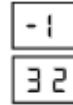

Seconds-mode


Press and hold "Start/STOP" button (C), and also press the "Power" button (A).

Seconds-mode display:


The time that was previously set will be stored, e.g.:

Note: max. 59 sec, speed range upto 6000 rpm.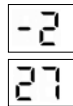

Setting the time

The time that was previously set will be stored, e.g.:
(Factory setting: 00)

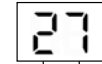Set the tens digit using button (E).

Set the ones digit using button (D).

Attaching the tube

Fit the tube carefully and ensure it is properly attached.

Setting the speed

Set the desired speed in the range of 300–6000 rpm using the infinitely adjustable rotary knob (B).

Starting

Press "Start/Stop" button (C), while the preset countdown is running, timer display will blink.

An audible signal will be given when the countdown is complete.

Cancelling the countdown

Press button "Start/Stop" (C).

Switching off

Press button "Power" (A).

Error message

It is possible for the media to become lodged between the cutters of the DT-tubes. The unit will automatically change the direction of rotation up to three times in an attempt to free the cutters. The unit will switch off automatically if this is not possible.

If the device becomes blocked on three occasions within a certain period of time the following error message will occur:

⇒ Display:

⇒ An audible signal will be given.


Corrective action: Switch off the device, check contents, switch on the device.

Accessories

- **ST-20** Tube with stirring device 20 ml, with pierceable membrane
- **ST-20-M-gamma** Tube with stirring device 20 ml, γ -sterilized, with pierceable membrane
- **ST-50** Tube with stirring device 50 ml
- **ST-50-M** Tube with stirring device 50 ml, with pierceable membrane
- **ST-50-M-gamma** Tube with stirring device 50 ml, γ -sterilized, with pierceable membrane
- **DT-20** Tube with dispersing device 20 ml, with pierceable membrane
- **DT-20 eco** Tube with dispersing device 20 ml, with PSU rotor/stator
- **DT-20-M-gamma** Tube with dispersing device 20 ml, γ -sterilized, with pierceable membrane
- **DT-50** Tube with dispersing device 50 ml
- **DT-50 eco** Tube with dispersing device 50 ml, with PSU rotor/stator
- **DT-50-M** Tube with dispersing device 50 ml, with pierceable membrane
- **DT-50-M-gamma** Tube with dispersing device 50 ml, γ -sterilized, with pierceable membrane
- **BMT-20-S** Tube for grinding with stainless steel balls 20 ml, with pierceable membrane
- **BMT-20-S-M-gamma** Tube for grinding with stainless steel balls, 20 ml, γ -sterilized, with pierceable membrane
- **BMT-50-S** Tube for grinding with stainless steel balls 50 ml
- **BMT-50-S-M** Tube for grinding with stainless steel balls 50 ml, with pierceable membrane
- **BMT-50-S-M-gamma** Tube for grinding with stainless steel balls 50 ml, γ -sterilized, with pierceable membrane
- **BMT-20-G** Tube for grinding with glass balls 20 ml, with pierceable membrane
- **BMT-50-G** Tube for grinding with glass balls 50 ml
- **BMT-50-G-M** Tube for grinding with glass balls 50 ml, with pierceable membrane
- **TC-20-M** Tube cover 20 ml, with pierceable membrane
- **TC-50** Tube cover 50 ml
- **TC-50-M** Tube cover 50 ml, with pierceable membrane
- **Glass balls**
- **Stainless steel balls**

Maintenance

The drive unit is maintenance-free but not immune from wear. The carbon brushes of the motor wear down over the time.

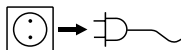For cleaning disconnect the mains plug.

Only use cleansing agents which have been recommended by **IKA®**.

Dirt

Dyes

Construction materials

Cosmetics

Foodstuffs

Fuels

Cleaning agent

isopropyl alcohol

water containing tenside/isopropyl alcohol

water containing tenside/isopropyl alcohol

water containing tenside

water containing tenside

For materials which are not listed, please request information from **IKA®** application support.

Wear protective gloves during cleaning the instruments.

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.

Do not allow moisture to get into the instrument when cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

Spare parts order:

When ordering spare parts, please give:

- machine type
- manufacturing number, see type plate
- item and designation of the spare part, see **www.ika.com**, spare parts diagram and spare parts list.

Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

For repair, please request the “**Decontamination Certificate**” form **IKA®** or use the download printout of it from **IKA®** website: **www.ika.com**.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

Technical data

Power supply unit

Input	VAC A Hz	100–240 0.8 50–60
Output	VDC Hz	24 24 W.LPS (limited power source) 50–60
Protection class		2 (double insulated) 

Tube Drive

Operating voltage	VDC mA	24 800
Power consumption, normal operation	W	20
Power consumption, standby operation	W	2
Power output	W	17
Drive		EC flat motor
Speed range: Mode 1 Mode 2	rpm	300 - 6000, infinitely adjustable up to 4000 up to 6000
Speed setting		rotating/pressing knob
Speed display		scale 0–9
Timer: mode 1 mode 2		1–29 min 1–59 sec
Display-timer		digital
Perm. ambient temperature	°C	+5 ... +40
Perm. rel. humidity	%	80
Protection to DIN 60529		IP 20
Contamination level		2
Overvoltage category		II
Operation at a terrestrial altitude level	m	max. 2000
Dimensions (W x D x H) without tube	mm	100 x 160 x 40
Weight	kg	0.75

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IKA®-Werke GmbH & Co.KG